

ScyTek
Laboratories

Instructions For Use
CEA-IFU

Rev. Date: March 4, 2009

Revision: 3

Page 1 of 1

P.O. Box 3286 - Logan, Utah 84323, U.S.A. - Tel. (800) 729-8350 – Tel. (435) 755-9848 - Fax (435) 755-0015 - www.scytek.com

Cresyl Echt Violet Solution (0.1%)

Description: Designed to show selective staining of Nissl granules. Nissl granules are purple, nuclei of neuroglia and endothelial cells are slightly bluer than Nissl granules.

Availability/Contents:

<u>Item #</u>	<u>Volume</u>
CEA125	125 ml
CEA500	500 ml
CEA999	1000 ml

Storage: 2-8° Centigrade

Procedure:

1. Deparaffinize sections if necessary and hydrate to distilled water.
2. Incubate slide in Cresyl Echt Violet (0.1%) for 2-5 minutes.
3. Rinse quickly in 1 change of distilled water.
4. Dehydrate quickly in 3 changes of absolute alcohol.
5. Clear, and mount in synthetic resin.

Storage: 2° C



8° C



ScyTek Laboratories, Inc.
205 South 600 West
Logan, UT 84321
U.S.A.



 EmergoEurope (31)(0) 70 345-8570
Molsstraat 15
2513 BH Hague, The Netherlands