



Instructions For Use TSS-IFU

Rev. Date: 10/29/03

Revision: 2

Page 1 of 1

P.O. Box 3286 - Logan, Utah 84323, U.S.A. - Tel. (800) 729-8350 - Fax (435) 755-0015 - www.scytek.com

Trypsin (Two Component Solution)

Description: Trypsin is a commonly used digestive enzyme for immunohistochemical procedures. This product is provided as a two component, ready-to-use solution.

Availability:

<u>Item #</u>	<u>Volume</u>
TSS155	30 ml Concentrate 125 ml Buffer
TSS310	60 ml Concentrate 250 ml Buffer

Storage: 2- 8 Centigrade

IMMUNOHISTOCHEMICAL Staining Protocol:

1. Combine trypsin concentrate (1 part), with buffer (3 parts) for a 0.125% trypsin solution. For stronger digestion, increase ratio to 1:2 or 1:1.
2. Incubate tissue section for 10 minutes at 37° Centigrade.
3. Rinse and proceed with immunostaining.